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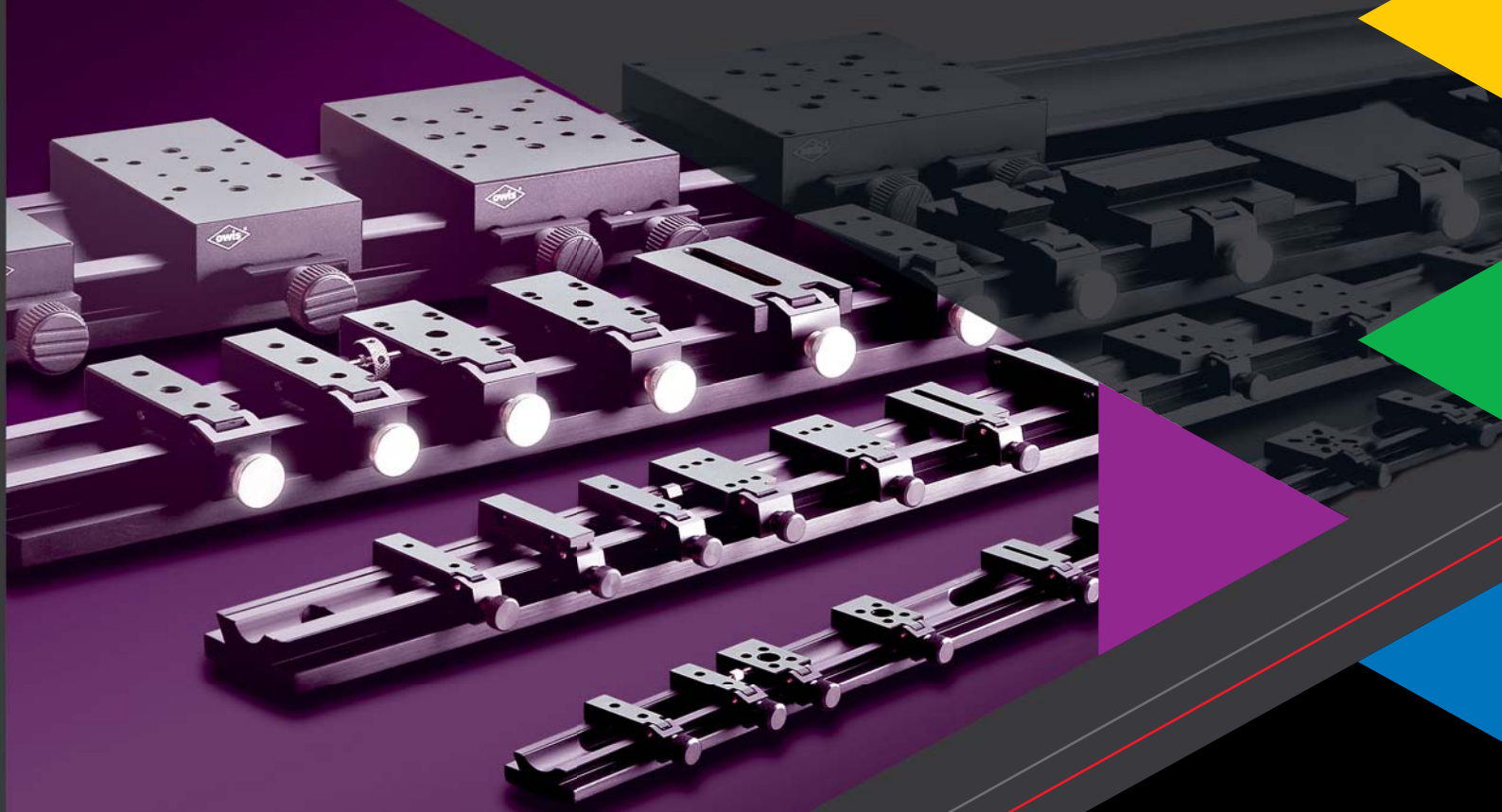
PRECISION IN PERFECTION

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OPTICAL BEAM HANDLING SYSTEMS

SYS 25
SYS 40
SYS 65
SYS 90

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SYS 25
SYS 40
SYS 65
SYS 90

*Hands-on training setups
Accessories*

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Content

Beam Handling Systems Accessories.....	ZB-S.....	S31
OWISet Hands-on training setups.....	PRAK.....	S29
Profiles, SYS 40.....	S 40-4.....	S9
Profiles, SYS 65.....	S 65-4.....	S17
Slides, SYS 25.....	RT 25.....	S3
Slides, SYS 40.....	RT 40.....	S11
Slides, SYS 65.....	RT 65.....	S19
Slides, SYS 90.....	RT 90.....	S25
System Rails with slotted holes, SYS 25.....	S 25 LL.....	S1
System Rails with slotted holes, SYS 40.....	S 40 LL.....	S7
System Rails with slotted holes, SYS 65.....	S 65 LL.....	S15
System Rails with slotted holes, SYS 90.....	S 90 LL.....	S23

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Symbol Legend:

In order to rapidly find the corresponding products, some symbols have been added to the product information.

-  for use with system SYS 25
-  for use with system SYS 40
-  for use with system SYS 65
-  for use with system SYS 90
-  optional non-magnetic versions
-  optional vacuum-prepared versions

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Type designation

PRAK.....OWISet Hands-on training setups.....	S29
RT 25SYS 25 Slides.....	S3
RT 40SYS 40 Slides	S11
RT 65SYS 65 Slides.....	S19
RT 90SYS 90 Slides	S25
S 25 LL.....SYS 25 System Rails with slotted holes.....	S1
S 40-4.....SYS 40 Profiles	S9
S 40 LLSYS 40 System Rails with slotted holes	S7
S 65-4.....SYS 65 Profiles	S17
S 65 LL.....SYS 65 System Rails with slotted holes.....	S15
S 90 LLSYS 90 System Rails with slotted holes	S23
ZB-S.....Beam Handling Systems Accessories.....	S31

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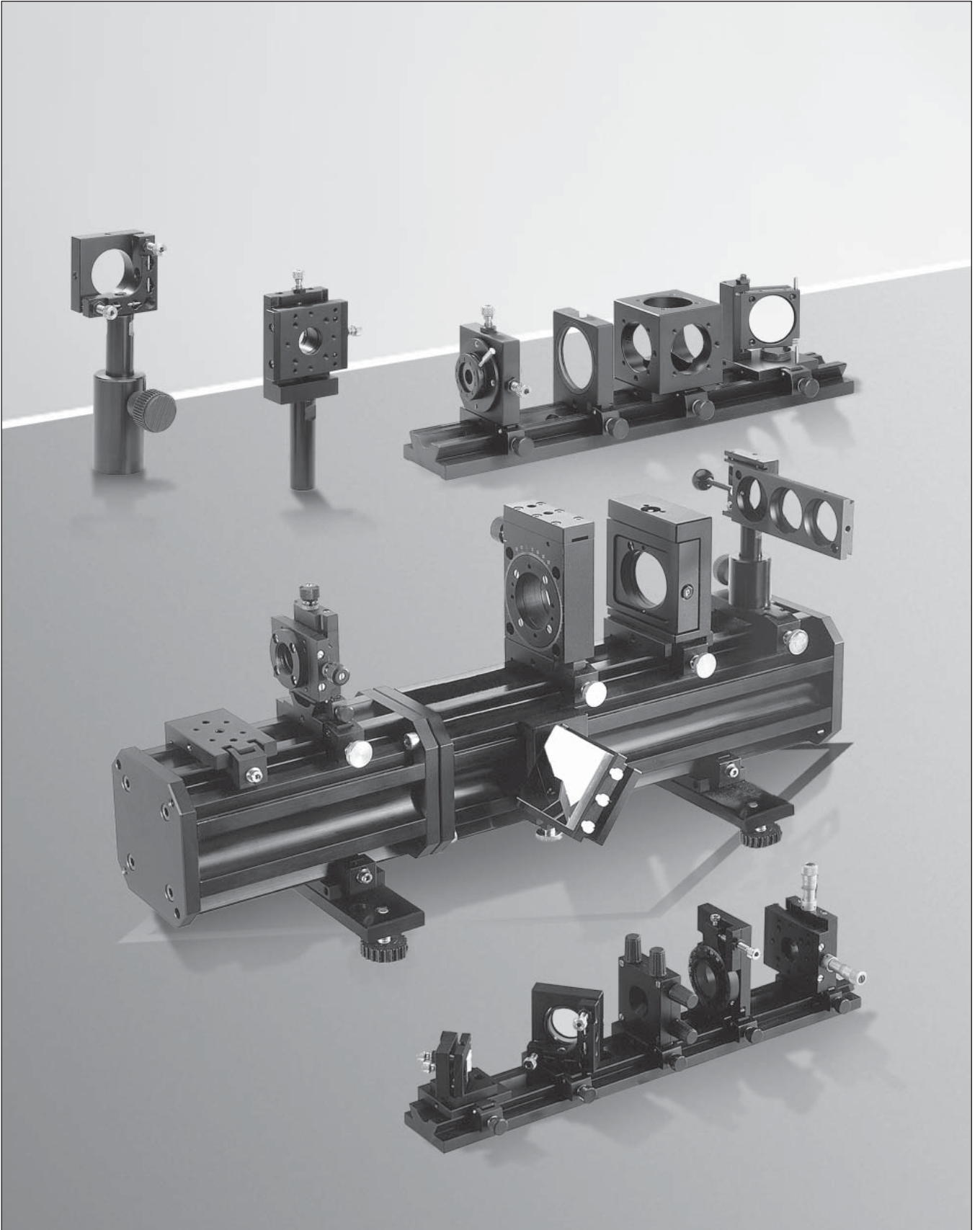
The OWIS® system installation kits allow the experimenter to provide experimental setups within shortest time. Furthermore, it is guaranteed that, when using the existing system components with a future series production, the same kits can be employed. This is particularly important for the industrial research. With the available system rails, the profiles and the corresponding riders, beam handling heights of 25 mm (SYS 25), 40 mm (SYS 40), 65 mm (SYS 65) and 105 mm (SYS 90) can be reached.

- rail/slide system allows rapid changing of components
- for linear, two- and three-dimensional setups
- versatile for laser and fibre-optic technology and classical and quantum optics
- for laboratory experiments and series productions
- SYS 25 components for use in SYS 40, SYS 40 components for use in SYS 65

Within the system, OWIS® tried to achieve the highest degree of compatibility among all system components.

A wide range of mounts and adapters make sure that all optical components can be properly mounted. Additionally these systems are compatible to systems from other manufacturers, who followed our standard.

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- andere Längen
- andere Eloxalfarben

Optionen

- other lengths
- other anodized colours

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The system rails S 25 LL are basic elements of the beam handling system SYS 25. They are available in standard lengths of 25 mm to 700 mm. The rails are made of deformation-resistant aluminium, precisely manufactured and reflection poor, black anodized.

The system rails, the slides and setup components of the SYS 25 system can be combined to stable and compact optical paths with an optical height of 25 mm.

Thanks to the slotted holes, the system rails S 25 LL can be mounted in almost any angle on the optical table. Thus, the variable handling required for optical setups is guaranteed.

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

Lieferung jeweils inkl. Schrauben ISO 7380 M6 x 10/screws ISO 7380 M6 x 10 are included in scope of delivery

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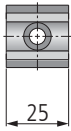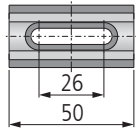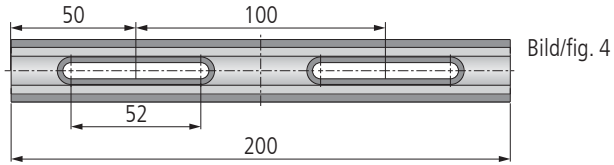
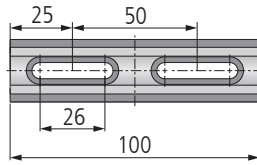
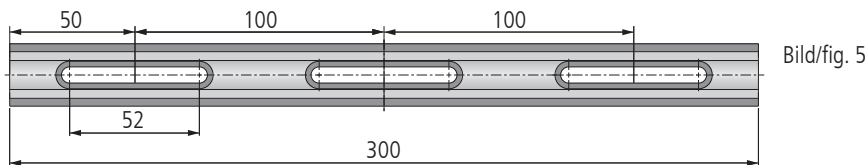
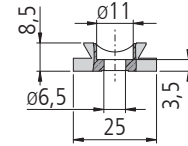
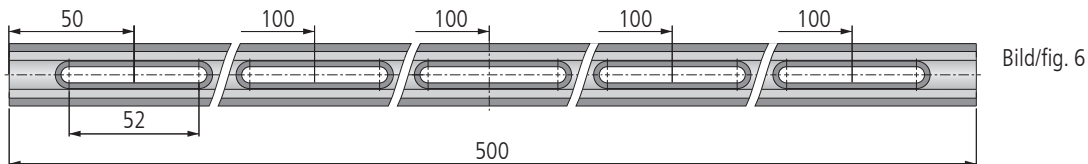
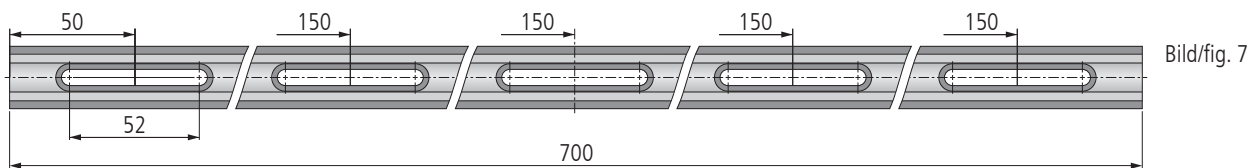
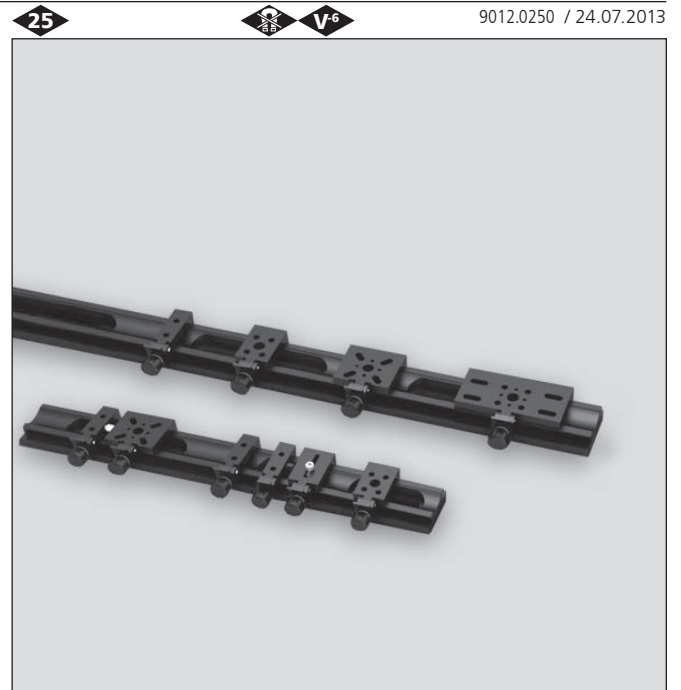
The RT 25 slides are used to mount assembling components on the system rail S 25 LL in order to provide the system height. The slides are equipped with the proven tilting clamping pieces. Thereby they can be placed easily also in existing setups on the system rails S 25 LL, too.



In order to fix the slides, a knurled or hexagon socket head cap screw is used. The form of the clamping pieces guarantees a high adhesion without constant deformation of the rail.

Slides with lateral threaded hole for the slide with fine adjustment on request.

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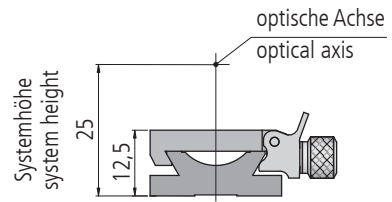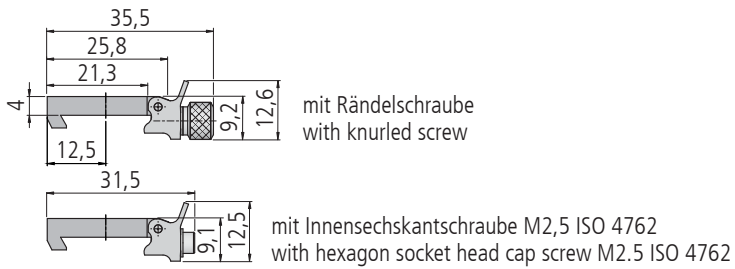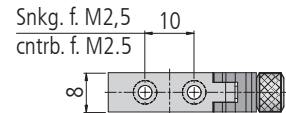
8 mm wide, for retaining mounting plates for lenses, filters etc.

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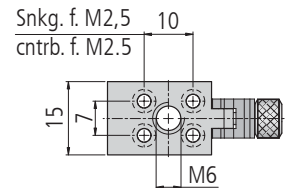
15 mm wide, for transmitting mounts, rotary mounts and mounting plates which need great stability and high precision.

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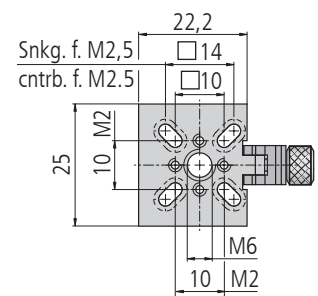
25 mm wide, for mounting mounting plates, optical mounts, for the other systems, too and for the VT 14.

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
incl. 4 slotted cheese head screws M2,5x5, 2 slotted cheese head screws M2x5

Typ/type	Bestell-Nr./part no.
RT 25-25-R	12.021.0025
RT 25-25-M2,5	12.022.0025

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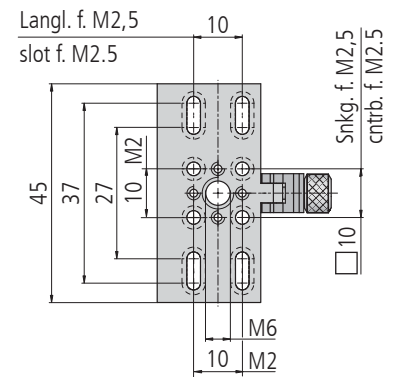45 mm wide, for mounting two single elements, mounting plates for example, or other large components.

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
incl. 4 slotted cheese head screws M2,5x5,
2 slotted cheese head screws M2x5

Typ/type	Bestell-Nr./part no.
RT 25-45-R	12.021.0045
RT 25-45-M2,5	12.022.0045

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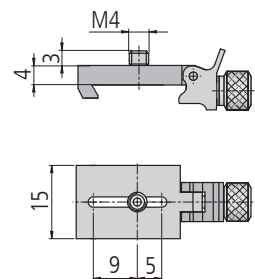15 mm wide, with slotted hole for mounting mirror and transmitting mounts in different angles.

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
incl. 1 slotted cheese head screws M2x4,
1 thread adapter M4/M2

Typ/type	Bestell-Nr./part no.
RT 25-15-R-LL	12.021.3015
RT 25-15-M2,5-LL	12.022.3015

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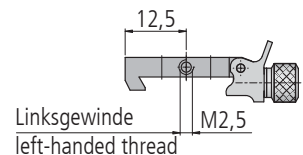
8 mm wide, with fine adjustment consisting of a slide with a left-handed frontal thread and a screw with right and left thread. Fits to slides with frontal M2.5 thread (on request). Max. adjustable range 6 mm.

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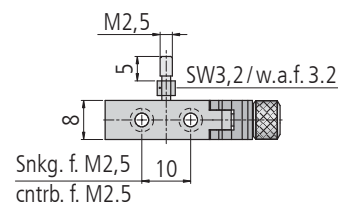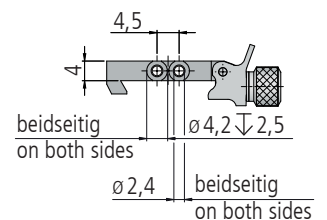
9 mm wide, with lateral counterbores for M2 to mount flip mirror holders KSH(M) 25 on S 25 LL. Two slides are included in scope of delivery.

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Optionen

- other lengths
- other anodized colours
- with tape scale

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The S 40 LL system rails are basic elements of the SYS 40 beam handling system. They are made of a special aluminium alloy, which ensures high dimensional stability. The guidance and assembly surfaces are machined with high precision. A sophisticated system concept allows precise and reproducible positioning of other SYS 40 components on the S 40 LL rails.

The system rails, RT 40 slides and other SYS 40 components, such as optic holders, form a system with a optical axis 40 mm above the base surface.

The slotted holes enable the mounting of the S 40 LL system rails on optical tables in virtually any angle to the thread grid.

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
Lieferung jeweils inkl. Schrauben ISO 4762 M6 x 12/screws ISO 4762 M6 x 12 are included in scope of delivery

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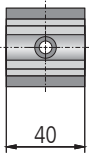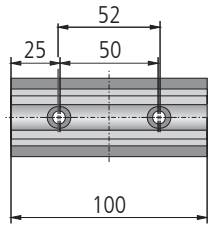
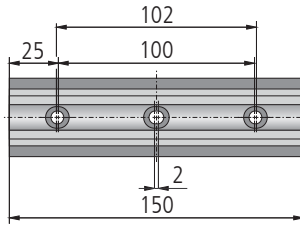
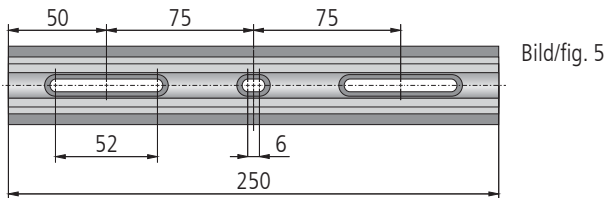
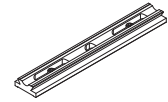
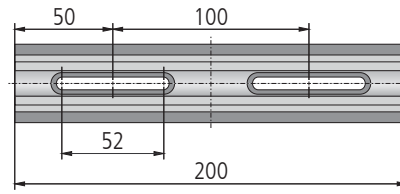
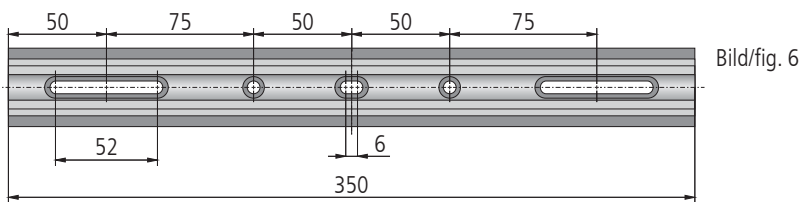
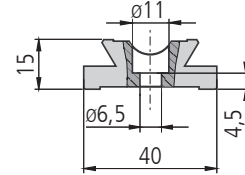
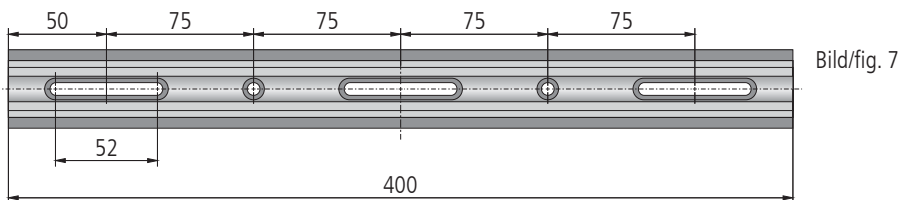
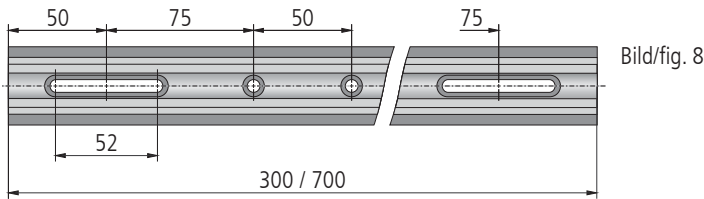
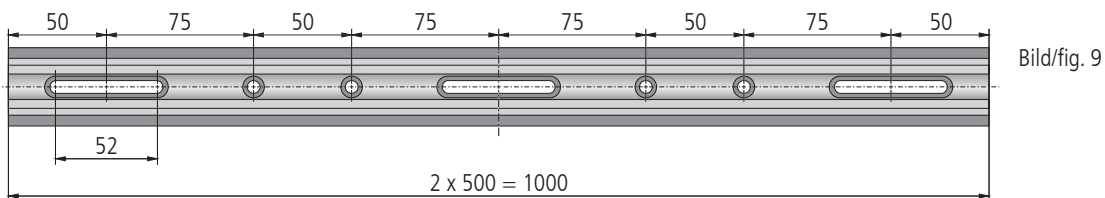
The 4-sided S 40-4 system profiles are the ideal choice for building 3D setups which are required to have a high dimensional stability. Their stiffness in torsion and bending even allows the building of cantilever constructions.

The four dovetail guides of the S 40-4 system profiles are fully compatible with those of the S 40 LL system rails. Thus, all SYS 40 equipment designed for the S 40 LL rails can be mounted on the S 40-4 system profiles as well.

The four guide rails of a profile are arranged in a way that allows four slides to be positioned around the profile independently of each other and without interference.

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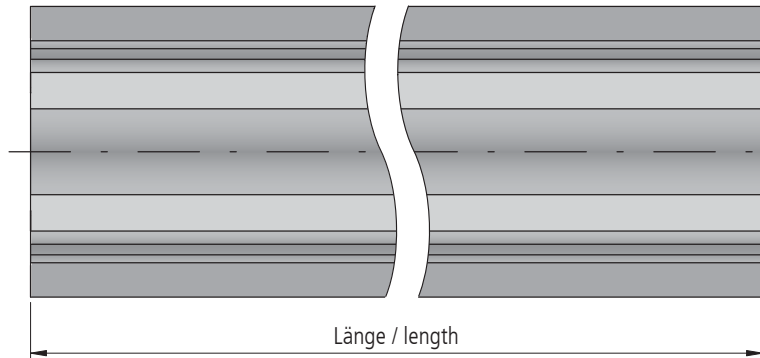
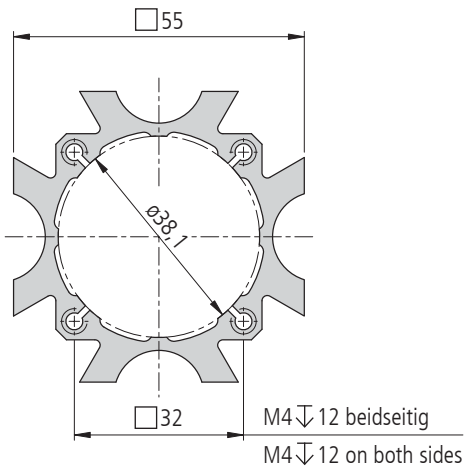
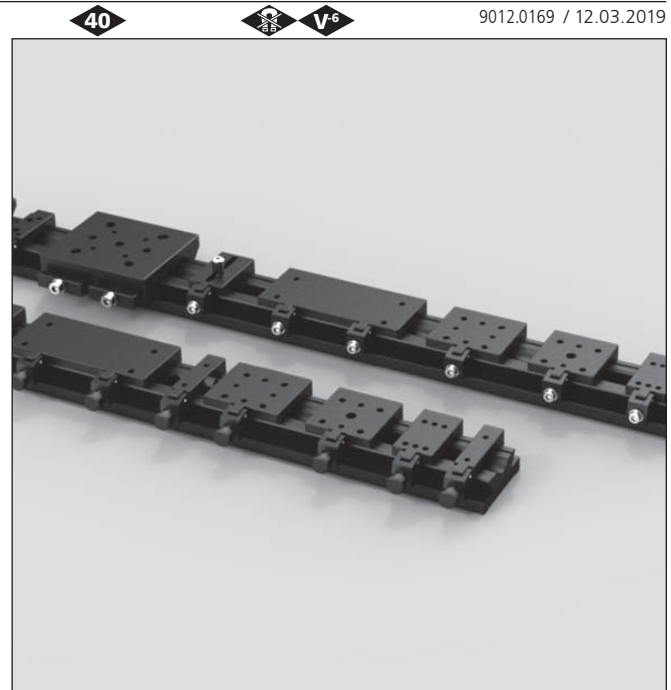
The RT 40 slides are equipped with tilting clamping pieces. The slides can be placed easily on the system rail S 40 LL or on the S 40-4 4-side profile and fixed with the screw. The structure of the clamping pieces guarantees a high adhesion without constant deformation of rail.

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

The RT 40 slides are available with knurled screw or with hexagon socket head cap screw ISO 4762.

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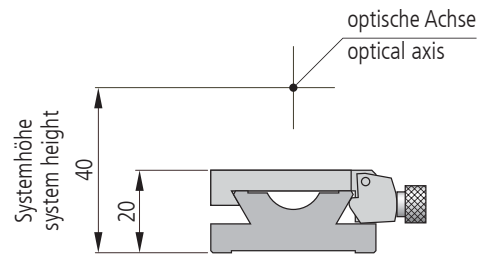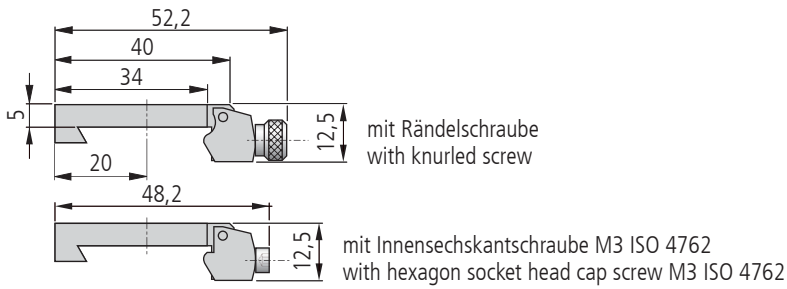
for retaining mounting plates for lenses,
filters etc.

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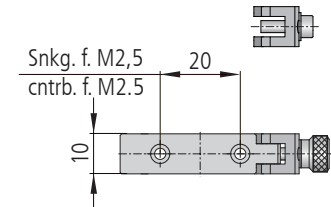
for transmitting mounts, rotary mounts and
mounting plates which need great stability and
high precision.

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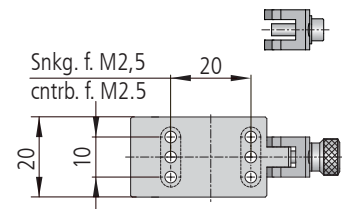
for mounting short pins or SL 12/16 columns
with M6 male threads, miniature XY stages
(horizontal), or translation stages.

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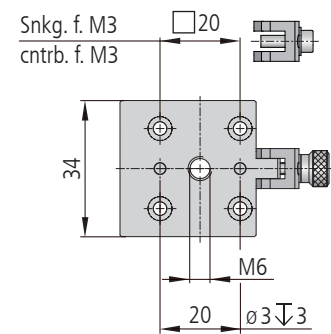
for mounting cubes, mounting plates,
miniature XY stages, optical mounts etc.
rotated 90° (e.g. on profile S 40-4).

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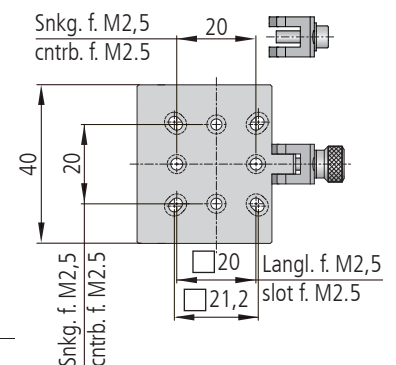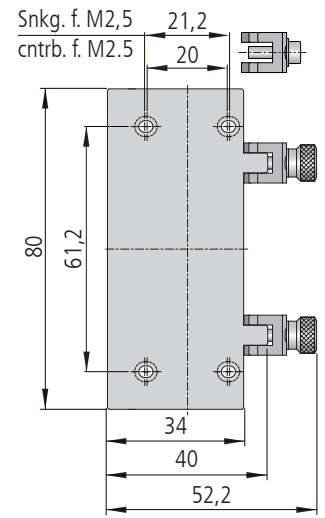
for mounting double cubes, other large components or two single elements, e.g. mounting plates.

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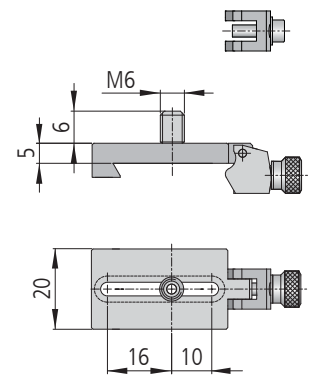
with slot for M3 screws for mounting mirror and transmitting adjustments. A grub screw with M6 external and M3 internal threads, is included.

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
incl. 1 hexagon socket head cap screw M3x8,
1 thread adapter M6/M3

Typ/type	Bestell-Nr./part no.
RT 40-20-R-LL	14.021.3020
RT 40-20-M3-LL	14.022.3020

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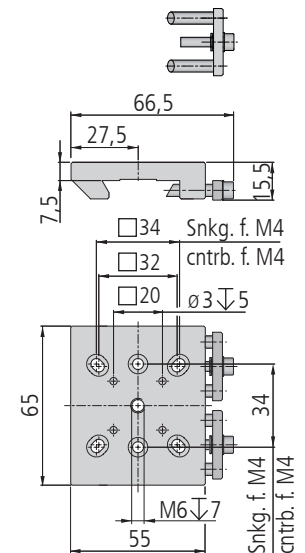has greater stability. It is used for building portals using S 40-4 profiles mounted on the front side. It also provides a stable mounting of system kit SYS 65 components if the system height is not really necessary.

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
incl. 4 hexagon socket head cap screws M4x6

Typ/type	Bestell-Nr./part no.
RT 40-4-65-M4	14.024.0065

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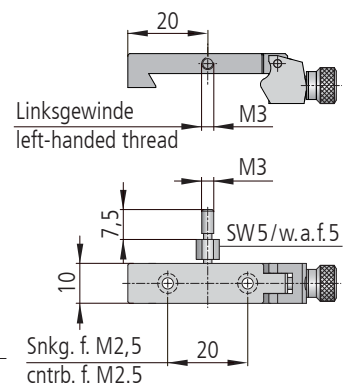fine adjustment consisting of a slide with a left-handed frontal thread and a screw with right and left thread. Fits to slides with frontal M3 thread (on request). Max. adjustable range 10 mm.

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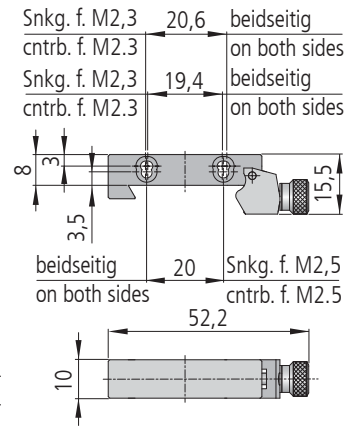
for horizontal mounting of WV 40 and to mount flip mirror holders KSH(M) 40 on S 40 LL. Two slides are included in scope of delivery.

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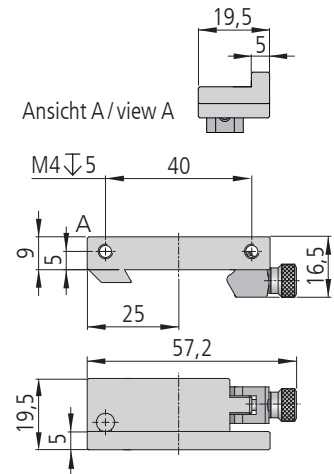
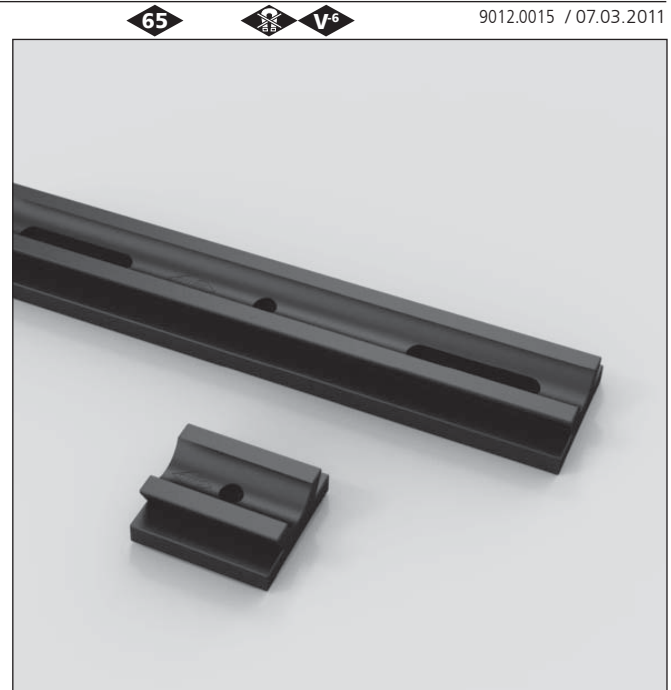
The S 65 LL system rails are basic elements of the SYS 65 beam handling system. They are made of a special aluminium alloy, which ensures high dimensional stability. The guidance and assembly surfaces are machined with high precision. A sophisticated system concept allows precise and reproducible positioning of other SYS 65 components on the S 65 LL rails.

The system rails, RT 65 slides and other SYS 65 components, such as optic holders, form a system with a optical axis 65 mm above the base surface.

The slotted holes enable the mounting of the S 65 LL system rails on optical tables in virtually any angle to the thread grid.

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
Lieferung jeweils inkl. Schrauben ISO 4762 M6x12/screws ISO 4762 M6x12 are included in scope of delivery

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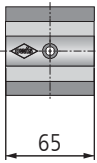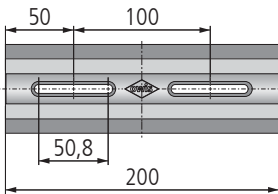
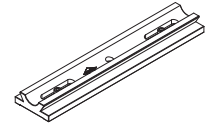
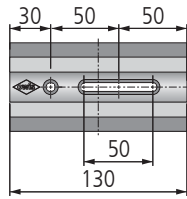
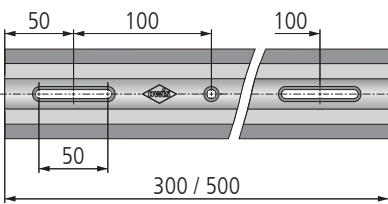
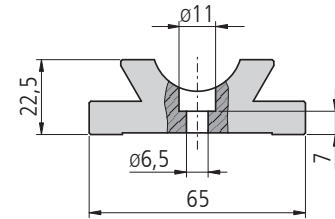
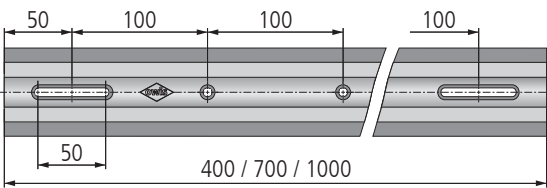
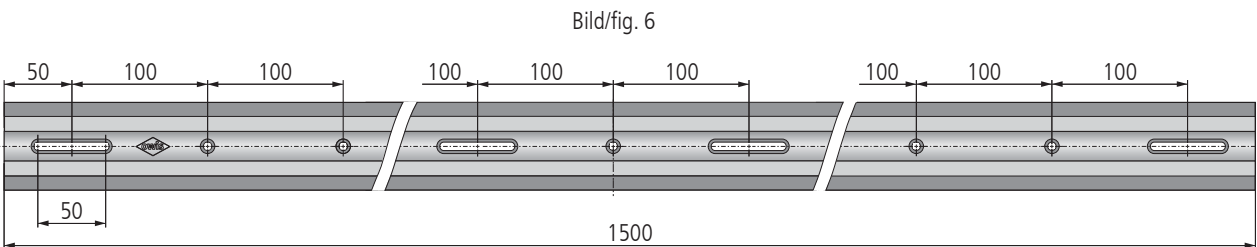
The 4-sided S 65-4 system profiles are the ideal choice for building 3D setups which are required to have a high dimensional stability. Their stiffness in torsion and bending even allows the building of cantilever constructions.

The four dovetail guides of the S 65-4 system profiles are fully compatible with those of the S 65 LL system rails. Thus, all SYS 65 equipment designed for the S 65 LL rails can be mounted on the S 65-4 system profiles as well.

The four guide rails of a profile are arranged in a way that allows four slides to be positioned around the profile independently of each other and without interference.

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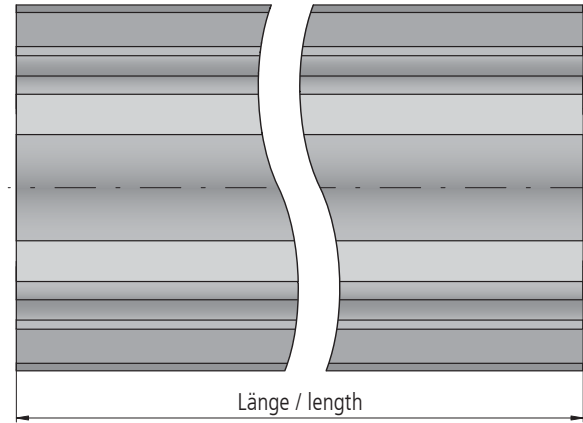
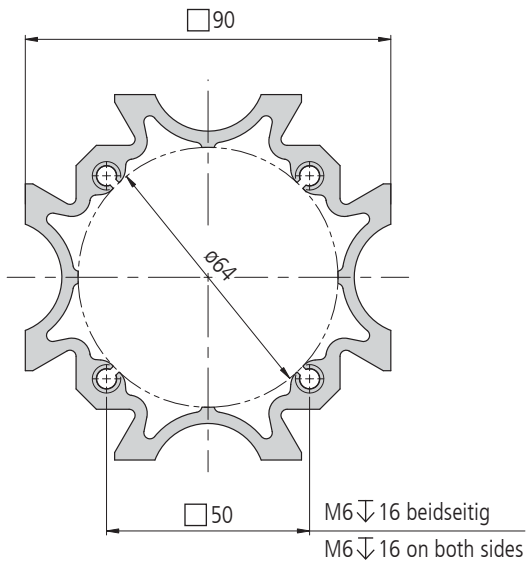
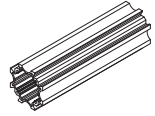
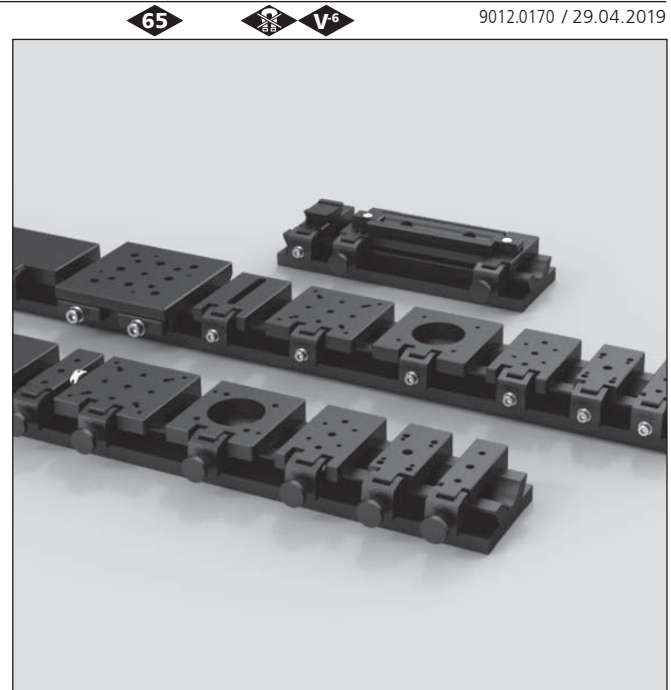
The RT 65 slides are equipped with tilting clamping pieces. The slides can be placed easily on the system rail S 65 LL or on the S 65-4 4-side profile and fixed with the screw. The structure of the clamping pieces guarantees a high adhesion without constant deformation of rail.

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

The slides RT 65 are available with knurled screw or with hexagon socket head cap screw ISO 4762.

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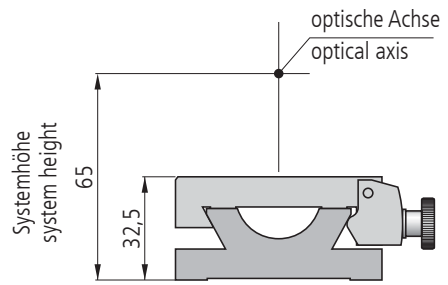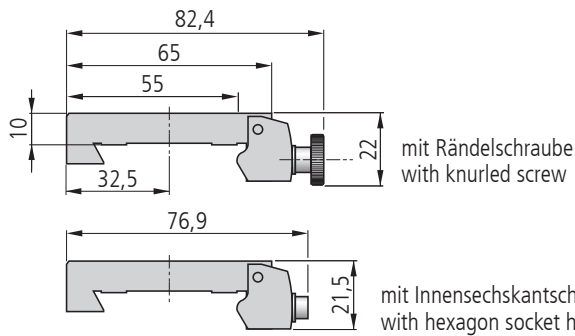
20 mm wide, for mounting carrier plates and other components with a 34 mm M4 grid.

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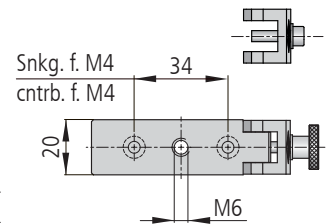
40 mm wide, for components with increased demands for stability and guidance, for components with 20 mm, M3 quadratic bore grids and/or with 34 mm M4 bore grids, as well as for columns with M6 base threads.

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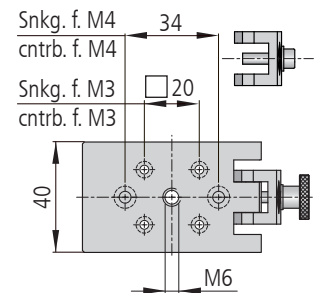
65 mm wide, additionally countersunk for M4 in a quadratic 34 mm and for M3 in a 38.9 mm grid and with M4 interior threads in a 50 mm grid.

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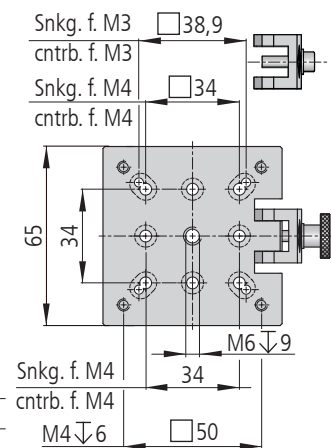
with slotted holes for M6 cylinder head bolt to mount mirror and optical mount at various angles.

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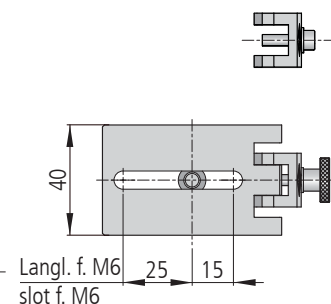
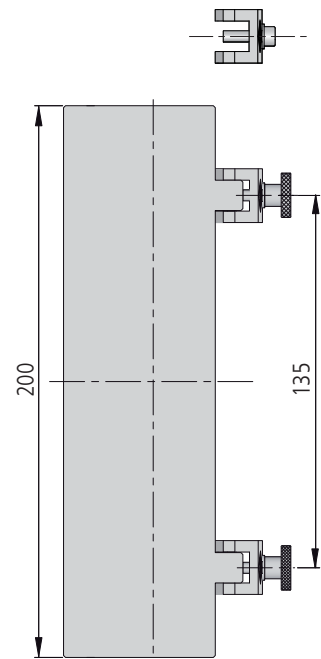
Experiment platform 65 for holding samples to be illuminated and examined with light. It can also hold laser cavities, optical cells, Q-switches, etc.. Two clamping devices ensure precise, reproducible positioning (also in axial direction in connection with a slide RT 65). The platform can thus be installed and removed easily during the experiment without maladjusting the rest of the setup.

Typ/type	Bestell-Nr./part no.
RT 65-200-R	16.021.0200
RT 65-200-M4	16.022.0200

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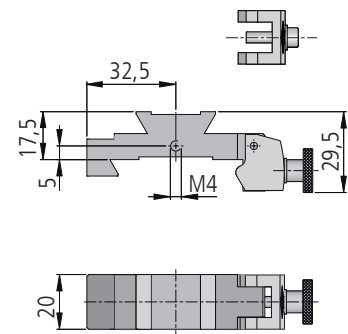
slide for SYS 65 with system rail S 40 LL, 20 mm, for adapting SYS 40 components to SYS 65. The optical height is therewith adjusted.

Typ/type	Bestell-Nr./part no.
RT 65-20-R-S40	16.091.4020
RT 65-20-M4-S40	16.092.4020

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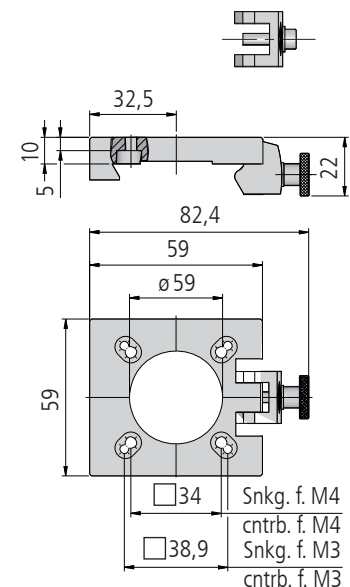Slide with central opening $\varnothing 35$ mm. For Cubes W 65 or for translation stages VT 65 or XY stages KT 65

Typ/type	Bestell-Nr./part no.
RT 65-59-R	16.021.1059
RT 65-59-M4	16.022.1059

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The S 90 LL system rails are basic elements of the SYS 90 beam handling system. They are made of a special aluminium alloy, which ensures high dimensional stability. The guidance and assembly surfaces are machined with high precision. A sophisticated system concept allows precise and reproducible positioning of other SYS 90 components on the S 90 LL rails.

The system rails, RT 90 slides and other SYS 90 components, such as optic holders, form a system with a optical axis 105 mm above the base surface.

The slotted holes enable the mounting of the S 90 LL system rails on optical tables in virtually any angle to the thread grid.

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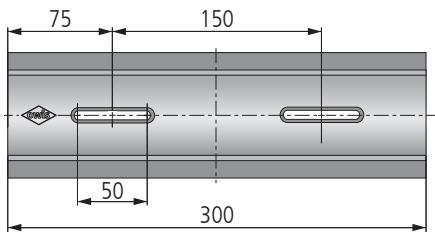
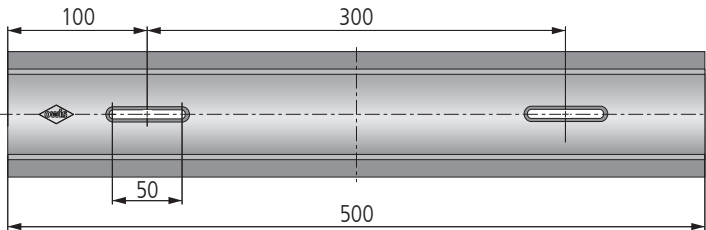
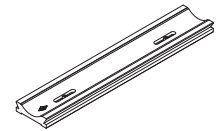
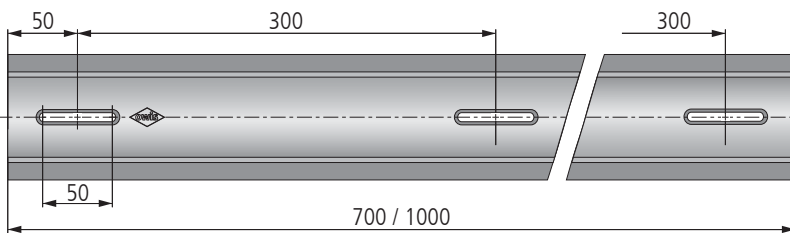
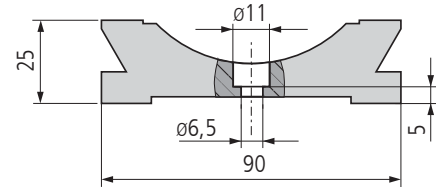
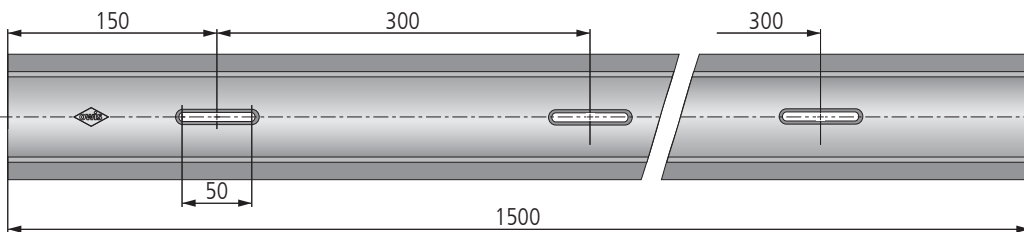
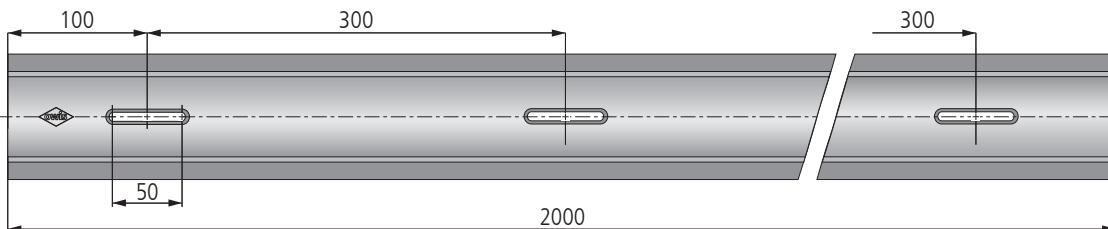
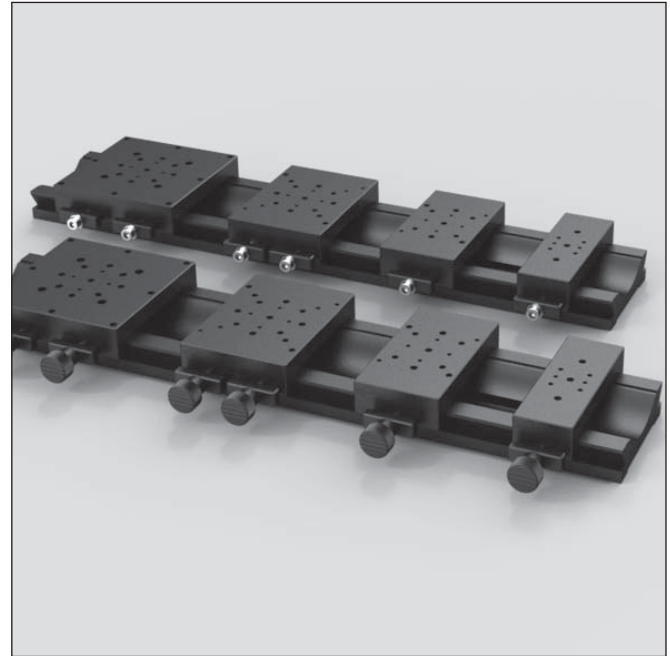
The RT 90 slides are used to realize setups with high beam heights. Also, they can carry large or heavy components.



To fix the RT 90 they are equipped with a clamping fork. These give a high stability to the slides. The structure of the clamping pieces guarantees a high adhesion without constant deformation of the rail.

The RT 90 slides are available with knurled screw or with hexagon socket head cap screw ISO 4762.

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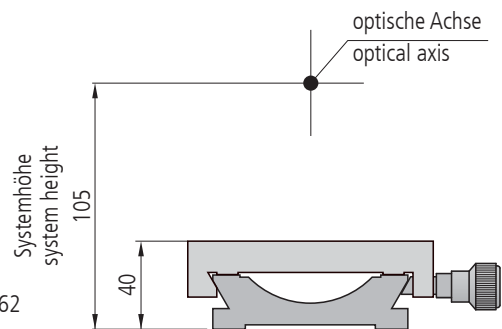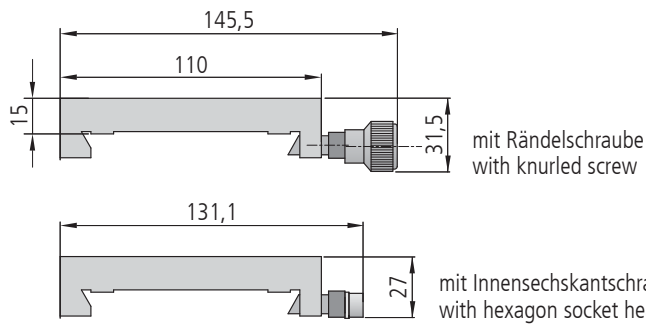
for fastening columns SL 12/16 with the highest possible stability. This lets you realize tightly spaced axial setups with large beam heights. Furthermore, the bore grids facilitate precise mounting of small measuring- and translation stages at right angles.


Typ/type	Bestell-Nr./part no.
RT90-40-R	19.021.0040
RT90-40-M6	19.022.0040

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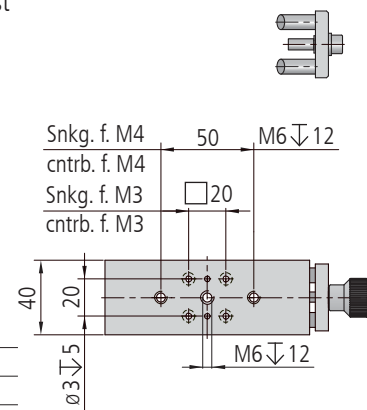for mounting large mirror- and transmitting mounts, columns with base plate, measuring- and translation stages, elevator stages, etc.


Typ/type	Bestell-Nr./part no.
RT90-60-R	19.021.0060
RT90-60-M6	19.022.0060

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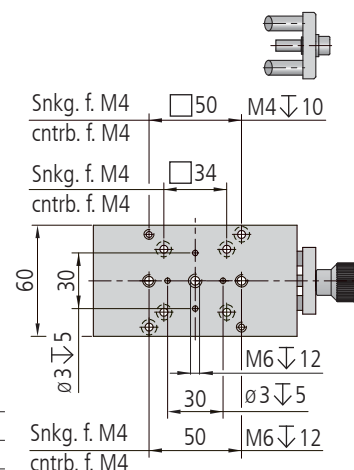
for an especially stable connection with a profile rail S90 LL through the use of two clamping forks. The bore grid includes five M6 threaded holes and an additional 70 mm M4 grid.

Typ/type	Bestell-Nr./part no.
RT90-80-R	19.021.0080
RT90-80-M6	19.022.0080

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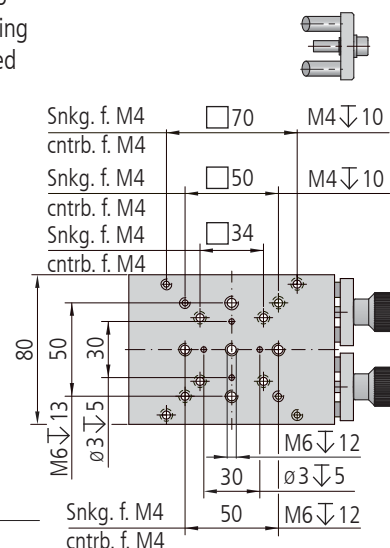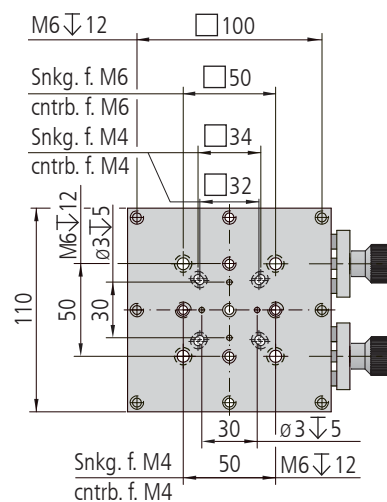for heavy loads and components with a M6 mounting grid, as well as for direct vertical mounting of S65-4 profiles.

Typ/type	Bestell-Nr./part no.
RT 90-110-R	19.021.0110
RT 90-110-M6	19.022.0110

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The OWISets PRAK for hands-on training setups are particularly suitable for use in institutions, universities and colleges. They provide all needed OWIS components for the hands-on experiments.

All sets are based on the beam handling system SYS 65. They can also be extended and combined with other OWIS components. If special setups are wanted /desired, the OWISets can also be individually configured or adapted.

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Adapter- and mounting plates are made of deformation-resistant aluminium. For protection, the surfaces are black anodized.

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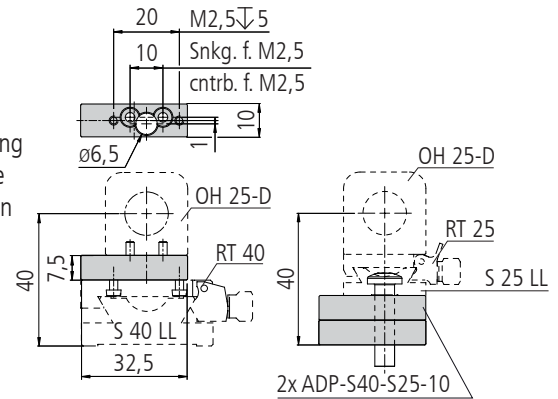
Adapter Plate SYS 40-SYS 25

are used to fix SYS 25 components on RT 40 slides as well as to adapt – by using two plates - the optical axes for a whole SYS 25 rail-slide-component combination to the beam height of the SYS 40.

incl. 2 screws ISO 1207 M2.5 x 6
incl. 2 screws ISO 4762 M2.5 x 8
incl. 1 screws ISO 7380 M6 x 25

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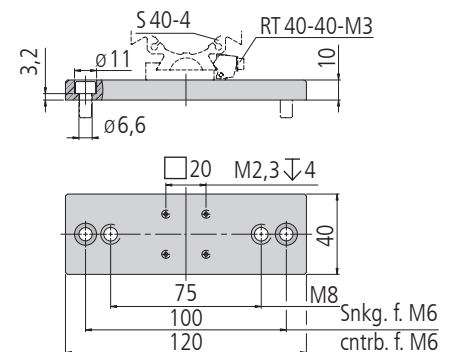
inkl. 2 Schrauben ISO 4762 M6 x 12
(nur MP 40-4-H)

Mounting Plate for Profile S 40-4, horizontal

is used to create optical benches or to mount frame constructions made of profiles on optical tables. Complete levelling feet based on the mounting plate are also available. They are connected on the bottom to the profile via slide RT 40, 40 mm wide.

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mounting plate for profile S 40-4, horizontal

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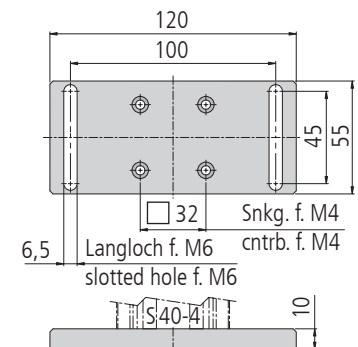
inkl. 4 Schrauben ISO 4762 M4 x 16
4 Schrauben ISO 4762 M6 x 16

Mounting Plate for Profile S 40-4, vertical

to mount vertically S 40-4 profiles on optical tables with 25 mm and 50 mm M6 grids. They are screwed down to the ends of the profiles via the four counterbores.

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mounting plate for profile S 40-4, vertical

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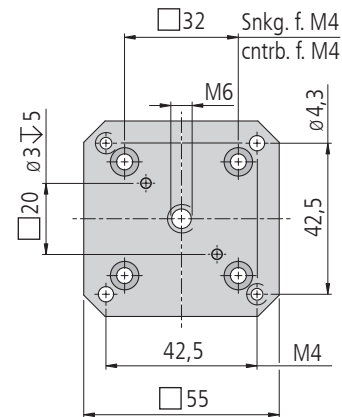
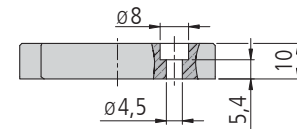
Connecting and End Plate for Profile S 40-4

is used to close off and stabilize the profiles, and is screwed down to the ends of the profiles. Furthermore, it makes it easy to connect two profiles together. To do this, the two opposing end plates are joined directly in 90° opposition to one another via the combined bore grid of two M4 threaded holes each and two holes for M4.

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
Bracket for Profile S 40-4

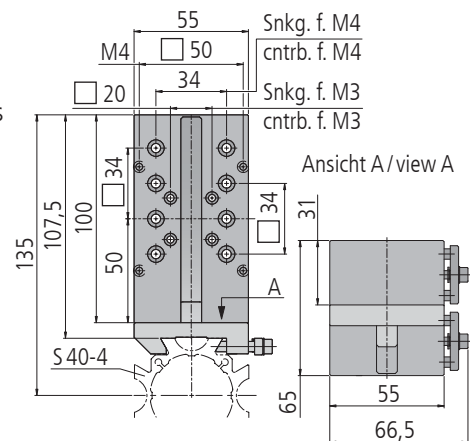enables to create horizontal mounting surfaces on vertically orientated profiles. It allows the mounting of measuring stages and translation stage combinations, prism stages, etc., with quadratic 20 mm bore grid for M3 and 34 mm for M4.

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
Rotary Base for Profile S 40-4 and S 65-4

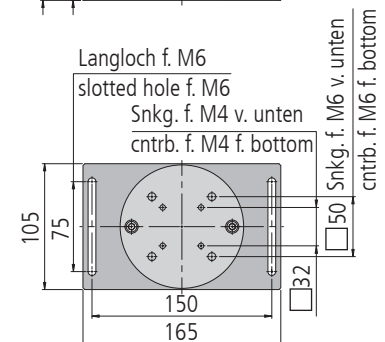make it simple to rotate columns by 90° around a definite pivot point on the profile axis without undermining stability.

inkl. 4 Schrauben ISO 4762 M4 x 16
4 Schrauben ISO 4762 M6 x 16
4 Schrauben ISO 4762 M6 x 25

incl. 4 screws ISO 4762 M4 x 16
4 screws ISO 4762 M6 x 16
4 screws ISO 4762 M6 x 25

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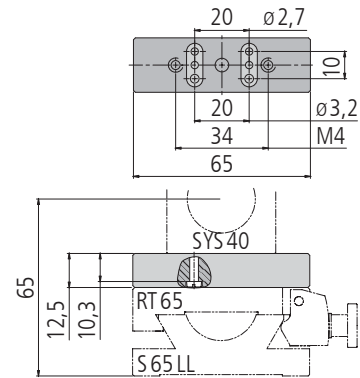
Adapter Plate SYS 65-SYS 40

to assemble SYS 40 components on RT 65 slides to adapt the optical axes.

incl. 2 screws ISO 1207 M2,3 x 14
2 screws ISO 1207 M2,5 x 14
2 screws ISO 1207 M3 x 16

	Typ/type	Bestell-Nr./part no.
Adapterplatte für SYS40-Komponenten in SYS 65	adapter plate for SYS 40 components in SYS 65	ADP-S65-S40-20
		14.094.0300

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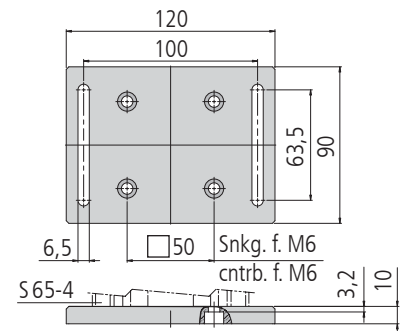
inkl. 8 Schrauben ISO 4762 M6 x 16

Mounting Plate for Profile S 65-4, vertical

to mount vertically profiles on optical tables with 25 mm and 50 mm M6 grids. It is screwed down to the ends of the profiles via the four counterbores.

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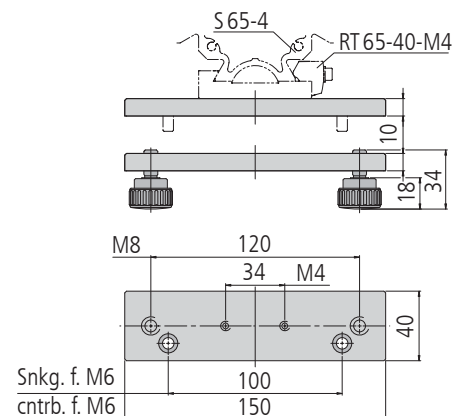
Mounting Plate / Levelling Foot, for Profile S 65-4, horizontal

is used to create optical benches or to mount frame constructions made of profiles on optical tables. Complete levelling feet based on the mounting plate are also available. They are connected on the bottom to the profile via slide RT 65-65 or slide RT 65-40.

incl. 2 screws ISO 4762 M6 x 12 (only MP 65-4-H)
2 adjusting screws M8 x 16 (only NF 65-4)


	Typ/type	Bestell-Nr./part no.
Montageplatte für Profil S 65-4, horizontal	mounting plate for profile S 65-4, horizontal	MP 65-4-H
		16.094.0100
Nivellierfuß für Profil S 65-4 mit Stellschrauben	levelling foot for profile S 65-4 with adjusting screws	NF 65-4
		16.094.0110

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Bracket for Profile S 65-4

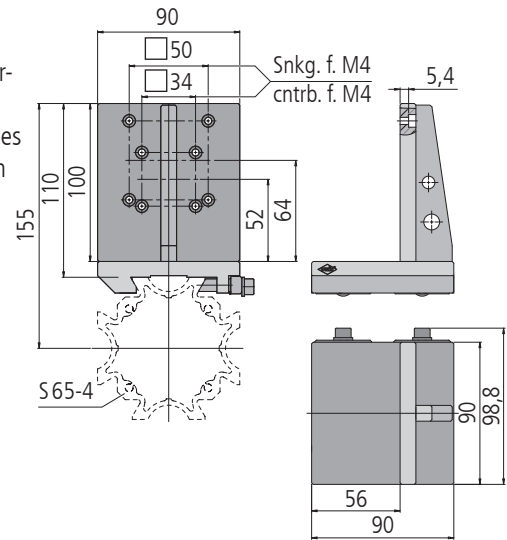enables to create horizontal mounting surfaces on vertically orientated profiles.
It allows the mounting of measuring stages and translation stage combinations, prism stages, etc., with quadratic bore grid 34 mm and 50 mm for M4.

inkl. 4 Schrauben ISO 4762 M4 x 12

incl. 4 screws ISO 4762 M4 x 12

		Typ/type	Bestell-Nr./part no.
Konsole für Profil S65-4	bracket for profile S65-4	KON-S65-4	16.094.0700