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Users are required to follow the safety and security advice.

Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

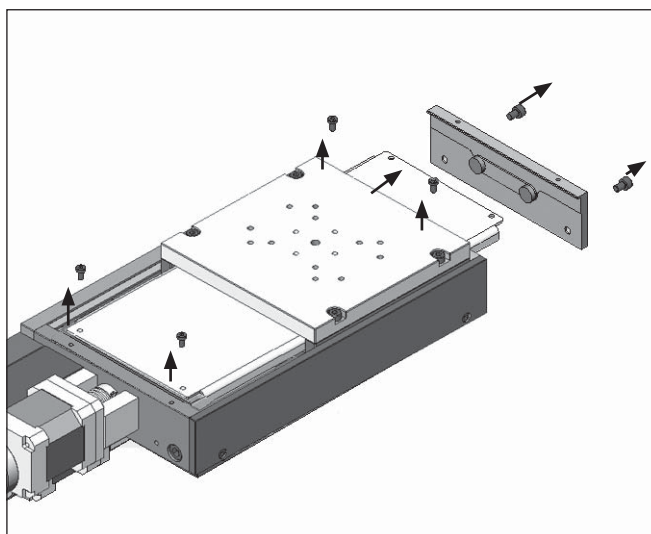
The **spindle** is lubricated by us with the **OWIS® SST.F11** grease.
Lubricating grease SST.F11 in a 5 ml applicator, part no. 90.999.0011

Relubrication should take place after about 500 operation hours, at least every 3 months.

The guides are maintenance free for 5 years.
Afterwards we suggest to send the product to OWIS® for an all-encompassing maintenance.

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Lubrication procedure

1. Drive the carriage to the end plate.
2. Turn off the control unit and make sure that it cannot be turned on accidentally.
3. Screw off the end plate.
4. Screw off the metal covering and slide it out backwards.
5. Remove the used grease or dust particles from the thread of the spindle resp. the ball screw with a dry, lint-free cloth. Cleansers such as benzene or petroleum are normally not necessary.
6. Apply evenly a thin film of lubricating grease SST.F11 onto the thread surface.
7. Insert the metal covering and screw it on.
8. Screw on the end plate.