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Users are required to follow the safety and security advice.

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Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

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The **worm gear** is lubricated by us with the **OWIS® SST.F8** grease. Lubricating grease SST.F8 in a 5ml applicator, part no. 90.999.0008

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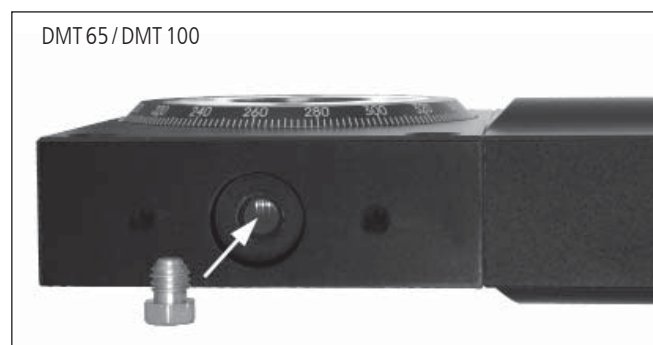
Relubrication should take place after about 500 operation hours, at least every 3 months.

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Lubrication procedure

1. Turn off the control unit and make sure that it cannot be turned on accidentally.
Attention! During the actual lubrication procedure the DMT must be halted, so that the applicator needle cannot be seized and pulled in by the worm.
2. Remove the lubricating nipple.
3. Insert the applicator needle into the open bore up to the worm and fill in about 0.1 ml grease SST.F8 (corresponds to ½ applicator scale-line).
4. Turn the plate another ¼ revolution and fill again about 0.1 ml grease SST.F8 into the opening. Continue this procedure until the rotary plate has accomplished a complete revolution.
5. Screw in the lubricating nipple.

Please note:

In the vacuum-prepared version no lubrication nipples are included. Therefore, steps 2 and 5 are omitted.

Furthermore, in vacuum-prepared versions, the grease SST.F5 must be used (instead of SST.F8)