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The JT 40T adjustable inserts enable optical elements to be set with high precision, secure against shock and vibration, and long-term stable. The moveable plate is supported on packets of disc springs, and held against the base plate by three screws.

Using an intermediate substrate of optical glass to absorb thermal stresses ensures a low-stress bond to eliminate distortion of the surface of the optical element.

This permits the use of adjustable inserts for beamsplitter cubes and plates, as well as front plane mirrors, or even for high-precision assemblies such as Mach-Zehnder or Michelson interferometers.

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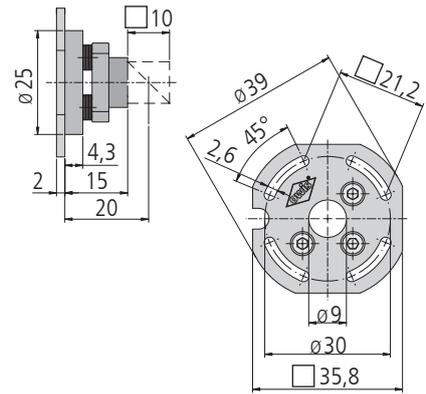
Adjustable Inserts for Beamsplitter Cubes and Prisms

with 10 mm or 15 mm side and optical glass substrate.

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
Adjustable Inserts for Beamsplitter Plates

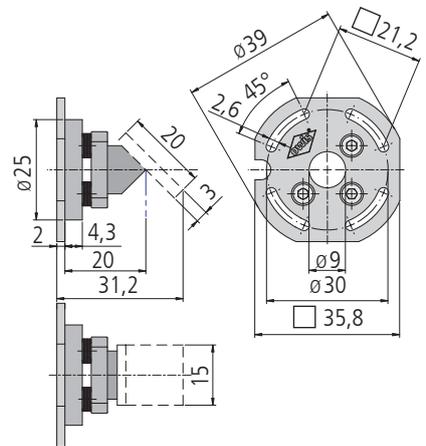with 15x20x3 mm and optical glass substrate.

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
Adjustable Inserts for Deflection Mirrors

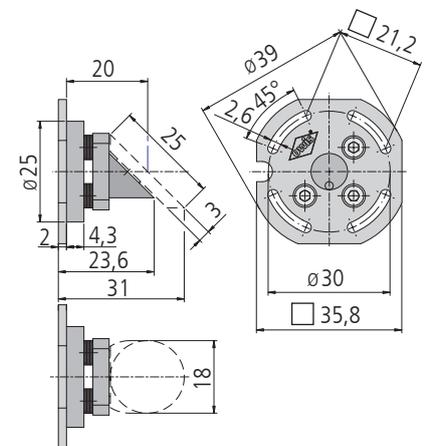with 18x25x3 mm for deflecting beams at 90°. Low-stress bonding to the optical glass substrate makes this suitable for high-quality laser mirrors. The optimum mirror thickness is 3 mm.

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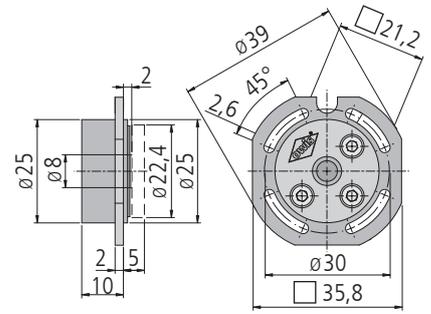
inkl. 4 Schrauben ISO 1207 M2,5x6

**Adjustable Inserts for Front Surface
plane and concave Mirrors**

with diameter 22.4 mm and optical glass substrate. The position of the mirror has been chosen so that a further adjustable insert, for example for a beamsplitter, can be installed in a cube W40.

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