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Users are required to follow the safety and security advice.

Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

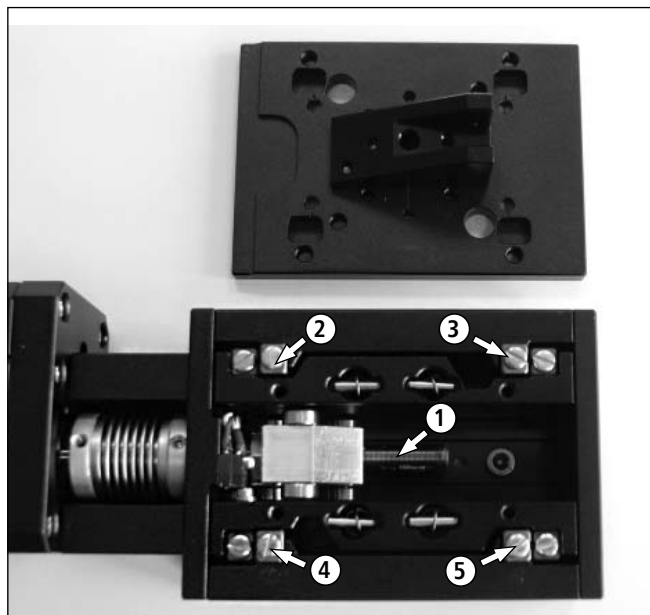
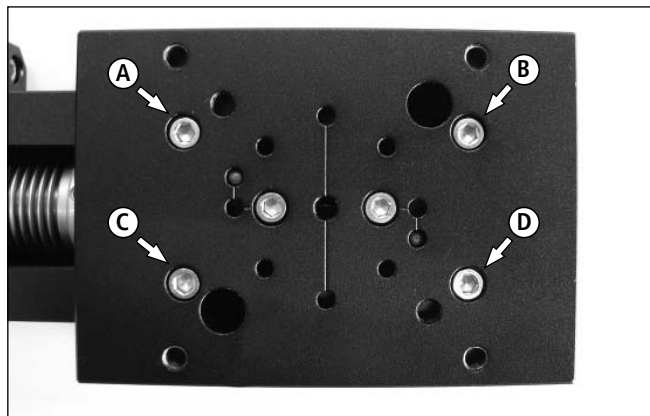
The **spindle** is lubricated by us with the **OWIS® SST.F11** grease.
Lubricating grease SST.F11 in a 5 ml applicator, part no. 90.999.0011

The **cross roller bearing guides** are lubricated by us with the **OWIS® SST.F1** grease.
Lubricating grease SST.F1 in a 5 ml applicator, part no. 90.999.0001

Relubrication should take place after about 500 operation hours, at least every 3 months.

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Lubrication procedure

1. Move the table plate into the lower end position.
2. Turn off the control unit and make sure that it cannot be turned on accidentally.
3. Screw off the table plate by removing the four screws (A-B).
4. Remove the used grease or dust particles from the threads of the fine-thread spindle (1) with a dry, lint-free cloth. Cleansers such as benzene or petroleum are normally not necessary.
5. Apply evenly a thin film of lubricating grease SST.F11 onto the thread surface (1).
6. Also apply a thin film of lubricating grease SST.F1 onto the rails (2), (3), (4) and (5) of the cross roller guides. For this purpose the guides have to be moved up and down manually.
7. Screw on the table plate. Consider the direction! It is important that the table plate is adjusted centrally and in parallel to all planes of the case.